POLYTRON® dispersing aggregates are made from stainless steel 1.4435 (316L), electric polished and can be used also for food processing.

The double bearing of the rotor shaft is a KINEMATICA standard, ensures quiet running, allows high rotor speeds up to 40'000 rpm and therewith it makes sure that extremely high tip speeds up to 50 m/s are reached for what POLYTRON® dispersing aggregates are known.

Heart of a POLYTRON® dispersing aggregate is always the rotor/stator dispersing head, called generator.

A speciality of KINEMATICA are technical sophisticated generator geometries for a wide range of homogenizing, dispersing, emulsifying, etc. applications.

Depending on the application and the product the most suitable dispersing aggregate can be chosen.

POLYTRON® dispersing aggregates of type G (with mechanical seal) are used for our REACTRON® RT 1 and RT 2 lab reactors. They ensure air tightness of the reactor vessel and allows working with vacuum.

The dispersing aggregates are driven by our standard lab **POLYTRON® dispersing drive units**.

Also your existing KINEMATICA drive units can be integrated in the REACTRON® systems.

The all-purpose drive systems PT 2100 and PT-K can be used for lab tasks from a few milliliters up to 2000 ml, respectively 3000 ml. They convince with easy handling, powerful motors and speeds up to 30000 rpm.

Together with the POLYTRON® dispersing aggregates, which are easily coupled to the drive unit without further tools, the product in the reactor vessel will be processed at the finest level in the micro scale range.

POLYMIX® stirrers, as a component of the REACTRON® system, ensure the macro mixing of the product in the reactor vessel. For our laboratory reactors our lab stirrer PX-SR 90 D has been chosen with the following features: large speed range, high torque, LED display, blocking and overheating protection.

Our REACTRON® systems can be **extended** to different degrees.

For example temperature control systems including thermostat / cryostat, temperature measuring and regulation or vacuum systems including vacuum pump and vacuum controller are available at KINEMATICA.

Call us. Our sales and applications department would be glad to assist you finding the best solution.

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Process Control Devices

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Finest Solutions ... worldwide!

A wide range of products with the finest solution for every application, matched by the superior quality of all components make KINEMATICA a world leader for Dispersing and Mixing technology.

With Keen Eyes ...

Development knows no rest at KINEMATICA. Our systems are constantly upgraded to reflect the state-of-the-art technology and the needs of the market. KINEMATICA has always had a keen eye for major advances, as proved by various patents. The rotor/stator principle for dispersing technology was conceived and developed at KINEMATICA and has become the world standard.

... For Your Needs

Applications for Dispersing/Homogenizing Technology are always on the increase. KINEMATICA continually creates new areas of processing applications based on innovations. We do not neglect the human factor as it is exemplified by the noise reduction of our high-speed systems.

Please, get in touch with us, our local appointed agent or your preferred dealer. We will be pleased to supply more details or arrange for a personal demonstration of our POLYTRON® and MEGATRON® systems.

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Represented by / Vertreten durch:

KINEMATICA AG
Dispersing- and Mixing Technology

Luzernerstrasse 147a
CH-6014 Littau / Lucerne
Switzerland

Tel.
Fax
e-Mail
web

+41-41-259 65 65
+41-41-259 65 75
laboratory@kinematica.ch
www.kinematica.ch

Printed in Switzerland (SIII) RT1-2 5/2003. Subject to technical modifications

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System REACTRON® RT 1 / RT 2 - Lab Reactors
Labor-Reaktoren

KINEMATICA AG
Dispersing- and Mixing Technology

REACTRON® Lab Reactor Systems / Labor-Reaktor Systeme

KINEMATICA is the leading manufacturer of homogenizing and mixing equipment based on our technology of rotor/stator dispersing principles.

Starting at smallest lab volumes of 0.1 ml up to production dimensions of 150 t/h we have the suitable equipment configuration on hand to solve your homogenizing, dispersing, emulsifying, suspending, wet milling and mixing task.

Our **POLYTRON®** and **MEGATRON®** homogenizing and mixing processors treats your product at the finest level in shortest time ("down sizing" and "micro mixing").

KINEMATICA is also a specialist for designing and manufacturing of complete plants. Our **REACTRON®** systems consist of processing tanks (single or multi-walled, standard design or complete customer and product specific), stirring system for macro mixing (e.g. anchor stirrer with drive unit), **POLYTRON® / MEGATRON®** homogenizing system for down sizing and micro mixing, piping, process control and further specific equipment in accordance with the customer, the product and the process.

REACTRON® systems are available for batch and semi-batch operation with volumes starting with 1 liter up to several 100 liters. They can be designed and manufactured according to customer and application specifications.

The **REACTRON® RT 1** and the **REACTRON® RT 2** systems are modular conceived lab reactors for optimizing and reproducing chemical, mixing, dispersing and homogenizing processes in lab scale.

Some application examples:

- Producing of cremes, lotions, emulsions
- Pharmaceutical or cosmetic products
- Suspending of solids in liquid media (e.g. liquid polymers)
- Dispersing of fine solids into a liquid or molten phase
- Suspending of additives and solid polymers in mineral oils
- Wet milling and dispersing of solids, fibers, stringy materials, tissues, cells into fluids and polymers
- Producing of chemical raw material and fine chemicals
- Producing of diary products and diet food
- Polymerisations
- Extracting of enzymes from bio masses
- *and further more*

Get in contact with our sales and applications departement to discuss your homogenizing, mixing, wet milling and complete plant tasks for lab, pilot plant and production. Together we will work out the *finest solution* to suit your application.

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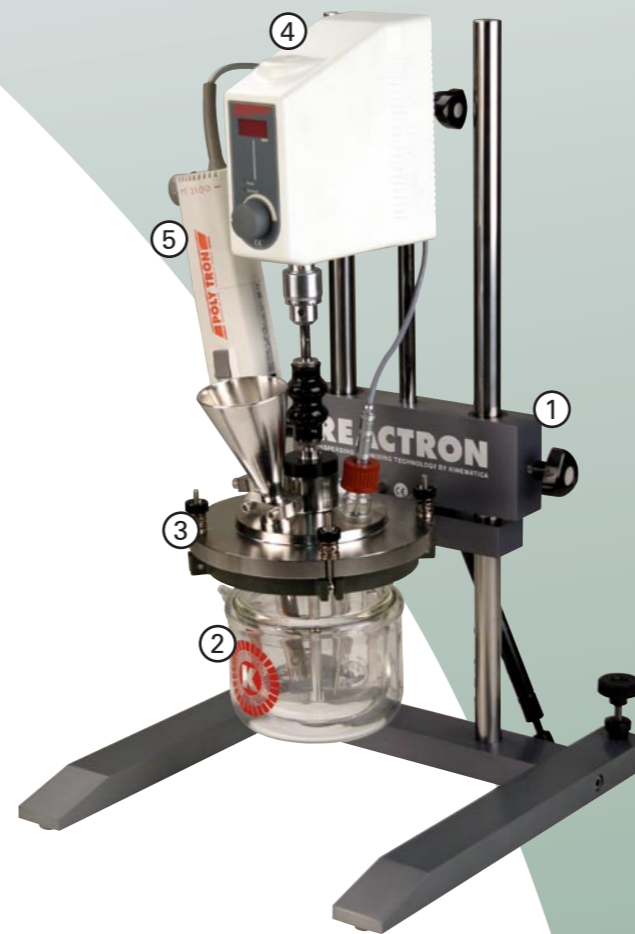
REACTRON® RT 150 (customer specific)
REACTRON® RT 150 (Kundenauslegung)

REACTRON®

DISPERSING AND MIXING TECHNOLOGY BY KINEMATICA

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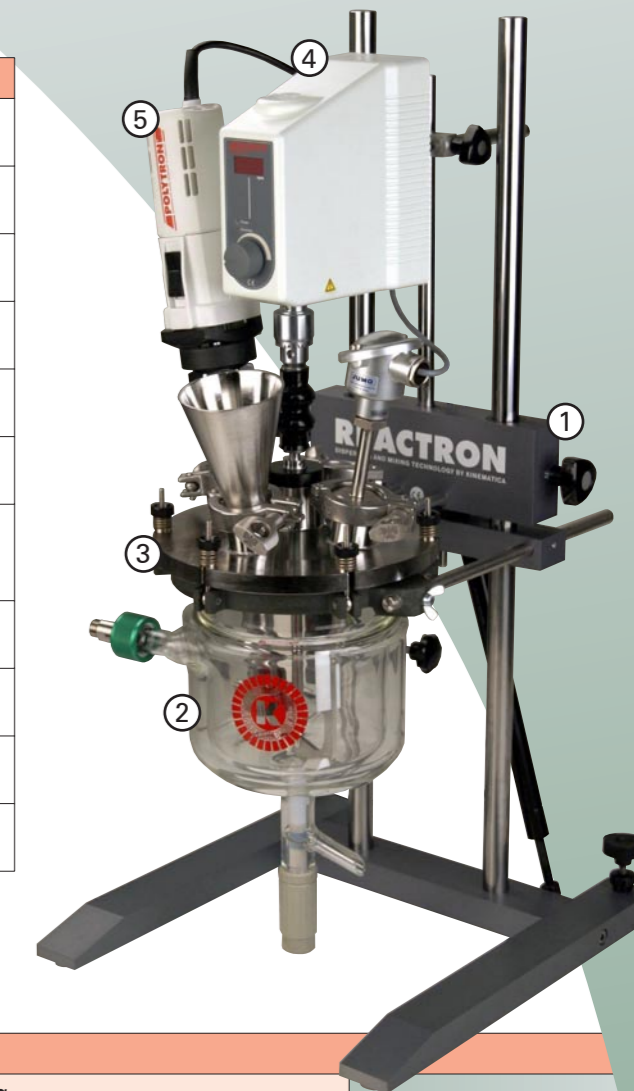
Dispersing Aggregate
PT-DA 2120/2G,
Anchor stirrer and
temperature sensor

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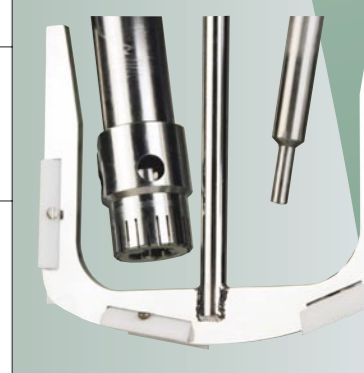
Dispersing Aggregate
PTA 30/4G,
Anchor stirrer and
temperature sensor

Dispergier-Aggregat
PTA 30/4G,
Ankerrührer und
Temperatursonde